

AppliClear - Water

Water bath disinfection

Product No. A7004

Description

AppliClear - Water is a concentrate for the prevention of growth of algae, moulds and has a germ-killing effect on microorganisms in water baths. It is bio-degradable and non-toxic.

Instructions:

The addition of 1 ml AppliClear - Water per 1 Liter water is sufficient to keep the water bath clean. This product is not suited for water baths in CO₂-incubators. AppliClear-Water is fully active, when the color indicator is blue. As soon as the color turns to colorless, refill with AppliClear-Water.

Storage: RT

AppliChem offers an alternative to AppliClear - Water:

Aquabator-Clean™ (100X)

Water bath disinfection

Product No. A9390

Description

Aquabator-Clean™ is intended for disinfecting various kinds of water baths from bacteria and fungi. It is recommended to use 10 ml Aquabator-Clean™ for each liter of water in the bath, and to repeat the procedure every 1 - 2 weeks. After 4 - 6 weeks, the bath should be emptied and refilled with water containing Aquabator-Clean™. No foam is produced when using the Aquabator-Clean™ with circulating water baths

The active ingredient in Aquabator-Clean™ is safe to humans and does not cause any irritating effects to the skin when used in the recommended concentration. It is biodegradable.

(Caution: Aquabator-Clean™ is not suited for use in CO₂-incubators!)

Storage: 2 - 8°C (protected from light)

Please note: If you need disinfectants for CO₂-incubators, we recommend the use of Incubator-Clean™ (A5230; spray for cleaning of the incubator) and Incuwater-Clean™ (A5219; concentrate for the water bath inside the incubator).